

BD Influx™ Cell Sorter

Filter Guide

Part #	Dyes	Application	# of Detectors	Filter(s)	Mirror(s)	Laser Blocking*
644708	355 nm, 100 mW + excitation optics					
644666	Hoechst	Cell cycle	1	460/50	N/A	400 LP
644667	Hoechst Blue Hoechst Red	Side population	2	460/50 670/30	550 DLP	400 LP
644668	Indo-1 bound Indo-1 free	Calcium flux	2	406/15 530/40	440 DLP	364 LP
650216	405 nm, 100 mW + excitation optics					
644669	Pacific Blue™ AmCyan	Cell surface markers	2	460/50 520/35	480 DLP	420 LP
646372	Pacific Blue™ Pacific Orange™	Cell Surface markers	2	460/50 550/50	480 DLP	420 LP
647490	Pacific Blue™ AmCyan Qdot® 655	Cell surface markers	3	460/50 520/35 650 LP	600 DLP	420 LP
646975	Pacific Blue™ AmCyan Qdot® 605 Qdot® 655/Qdot® 705	Cell surface markers	4	460/50 520/35 605/40 650 LP	480 DLP 575 DLP 645 DLP	420 LP
650198	BD Horizon™ V450 BD Horizon V500 Qdot® 605 Qdot® 655 Qdot® 705 Qdot® 800	Cell surface markers	6	460/50 520/35 610/20 660/20 710/50 750 LP	480 DLP 550 DLP 645 DLP 685 DLP 740 DLP	420 LP
644710	457 nm, 300 mW + excitation optics					
644670	CFP Chromomycin	Fluorescent protein, chromosomes	2	480/40 550/50	500 DLP	458 LP
644671	CFP Primary chlorophyll	Fluorescent protein, marine biology	2	480/40 650 LP	610 DLP	458 LP
646356	488 nm, 200 mW + excitation optics					
644672	FITC PE	Cell surface markers	2	530/40 580/30	550 DLP	505 LP
644673	GFP YFP	Fluorescent proteins	2	513/17 542/27	527 DLP	505 LP
644674	FITC/GFP Chlorophyll	Fluorescent proteins, marine biology	2	530/40 692/40	600 DLP	505 LP
644675	FITC PI	Cell surface markers, live/dead discrimination, cell cycle	2	530/40 610/20	550 DLP	505 LP
644676	FITC PerCP PerCP-Cy™5.5	Cell surface markers	3	530/40 670/30 710/50	550 DLP 685 DLP	505 LP
644677	FITC PE PE-Texas Red®/propidium iodide PE-Cy™5	Cell surface markers, live/dead discrimination, cell cycle	4	530/40 580/30 610/20 670/30	550 DLP 600 DLP 645 DLP	505 LP
644678	FITC PE PE-Cy5 PE-Cy™7	Cell surface markers	4	530/40 585/29 670/30 750 LP	550 DLP 610 DLP 710 DLP	505 LP
644679	FITC/SYBR® Green PE PE-Texas Red®/propidium iodide Chlorophyll	Cell surface markers, live/dead discrimination, cell cycle, marine biology, microbiology	4	530/40 580/30 610/20 692/40	550 DLP 600 DLP 645 DLP	505 LP
644680	FITC/SYBR® Green SYBR® Gold PE Chlorophyll	Cell surface markers, marine biology, microbiology	4	520/15 542/27 585/40 692/40	527 DLP 560 DLP 610 DLP	505 LP

DLP = Dichroic longpass LP = Longpass

* Depending on the combination of laser lines in an instrument configuration, more rigorous laser blocking might be necessary. BD Biosciences will add high-quality laser line blocking filters when necessary.



BD Influx™ Cell Sorter Filter Guide

Part #	Dyes	Application	# of Detectors	Filter(s)	Mirror(s)	Laser Blocking*
646356 488 nm, 200 mW + excitation optics						
644681	FITC	Cell surface markers, live/dead discrimination, cell cycle	5	530/40	550 DLP	505 LP
	PE			580/30	600 DLP	
	PE-Texas Red®/propidium iodide			610/20	645 DLP	
	PE-Cy5.5			692/40	740 DLP	
	PE-Cy7			750 LP		
644712 532 nm, 150 mW + excitation optics						
644682	Primary PE	Cell surface markers	2	585/40	645 DLP	532 LP
	PE-Cy5			670/30		
644683	PE	Cell surface markers	3	580/30	600 DLP	532 LP
	PE-Cy5			660/40	680 DLP	
	PE-Cy5.5			710/50		
644684	PE	Cell surface markers, live/dead discrimination, cell cycle	4	585/29	600 DLP	532 LP
	PE-Texas Red®/propidium iodide			610/20	645 DLP	
	PE-Cy5			670/30	700 DLP	
	PE-Cy7			750 LP		
644689	PE	Cell surface markers	4	593/40	645 DLP	570 LP
	PE-Cy5			670/30	685 DLP	
	PE-Cy5.5			710/50	740 DLP	
	PE-Cy7			750 LP		
644713 561 nm, 75 mW + excitation optics						
650217 561 nm, 150 mW + excitation optics						
644685	DsRed	Fluorescent proteins	2	585/29	600 DLP	570 LP
	mCherry			615/24		
644686	PE	Cell surface markers	2	593/40	645 DLP	570 LP
	PE-Cy5			670/30		
644687	PE	Cell surface markers	3	593/40	645 DLP	570 LP
	PE-Cy5			670/30	700 DLP	
	PE-Cy7			750 LP		
644688	PE	Cell surface markers	4	585/29	600 DLP	570 LP
	PE-Texas Red®			610/20	645 DLP	
	PE-Cy5			670/30	700 DLP	
	PE-Cy7			750 LP		
644689	PE	Cell surface markers	4	593/40	645 DLP	570 LP
	PE-Cy5			670/30	685 DLP	
	PE-Cy5.5			710/50	740 DLP	
	PE-Cy7			750 LP		
648728	PE	Cell surface markers	5	585/29	600 DLP	561 LP
	PE-Texas Red®			610/20	645 DLP	
	PE-Cy5			670/30	685 DLP	
	PE-Cy5.5			710/30	740 DLP	
	PE-Cy7			750 LP		
650218 640 nm, 120 mW + excitation optics						
644690	APC	Cell surface markers	2	670/30	700 DLP	650 LP
	APC-Cy7			750 LP		
644691	Chlorophyll/APC	Cell surface markers, marine biology	2	692/40	740 DLP	650 LP
	APC-Cy7			750 LP		
646371	APC	Cell surface markers	3	670/30	700 DLP	650 LP
	Alexa Fluor® 700			720/40	740 DLP	
	APC-Cy7			750 LP		

DLP = Dichroic longpass LP = Longpass

* Depending on the combination of laser lines in an instrument configuration, more rigorous laser blocking might be necessary. BD Biosciences will add high-quality laser line blocking filters when necessary.



Class I (1) laser product.

For Research Use Only.

Cy™ is a trademark of Amersham Biosciences Corp. Cy dyes are subject to proprietary rights of Amersham Biosciences Corp and Carnegie Mellon University and are made and sold under license from Amersham Biosciences Corp only for research and in vitro diagnostic use. Any other use requires a commercial sublicense from Amersham Biosciences Corp, 800 Centennial Avenue, Piscataway, NJ 08855-1327, USA.

Pacific Blue and Pacific Orange are trademarks and Alexa Fluor, Texas Red, and SYBR are registered trademarks of Molecular Probes, Inc.

BD, BD Logo and all other trademarks are property of Becton, Dickinson and Company. © 2011 BD

23-10507-04

BD Biosciences
2350 Qume Drive
San Jose, CA 95131
US Orders: 877.232.8995
bdbiosciences.com